

Jamieson Laboratories Relies on Impulse®/P Size Grader from Symetix®

SYMETIX®

As Canada's oldest and largest manufacturer of natural vitamins and minerals, Jamieson Laboratories is a market leader that has earned that position by consistently providing innovative products of the highest quality, purity, and safety. To help maintain their high standards, Jamieson recently installed an Impulse®/P high-capacity vibratory size grader from Symetix®. Jamieson's Impulse/P automates the removal of broken tablets and fragments prior to packaging to improve product quality while reducing labor costs.

"Prior to Impulse/P, we were manually sorting tablets for breakage. We searched for an automated solution that would improve product quality and alleviate unnecessary labor at the same time," said Andy Holwell, Director of Operations at Jamieson. "What used to take a week to manually sort on an inspection belt can now be accomplished in an eight-hour shift, including clean up."

Designed for solid-dose pharmaceutical and nutraceutical manufacturers and contract packers, the Impulse/P vibratory size grader features a sanitary stainless steel bed and removable decks

with holes that are perfectly sized for each product. Gentle electromagnetic power is combined with a high-throughput design to mechanically sort softgels, (system does not process two piece capsules) and tablets.

Jamieson selected Symetix's Impulse/P for tablets, a 48-inch (122-cm) wide system that size grades up to 1,000,000 tablets per hour, while their subsidiary, International Nutrient Technologies, installed the 18-inch (46-cm) wide Impulse/P for softgels, which size grades up to 500,000 softgels per hour. Jamieson's two-deck tablet system removes broken tablets and fragments, and the three-deck softgel system at International Nutrient Technologies removes twins, clusters, slugs, under-diameter capsules, and over-diameter softgels.

"Our goal was to achieve performance results that matched the sales pitch. Symetix told us that Impulse/P could remove 99 percent of broken tablets and fragments. We're getting 99.8 percent removal. That level of performance can't be accomplished with manual inspection, especially at those high speeds," said Adriana Bodica, Process Engineer. "The upper





deck removes broken tablets and fragments, and the lower deck recovers usable tablets, which helps keep nuisance rejects – the loss of good tablets – to a very low and acceptable volume.”

The decks are removable – each tablet length to be graded has a corresponding set of dedicated decks, which assures repeatable grading. Jamieson has three sets of decks to handle their range of products. “When we were validating Impulse/P, we did dozens of tests back-to-back to be sure the performance was consistent from one run to another. It is very repeatable,” noted Bodica.

Compared to a traditional diverging roller size grader, which introduces subjectivity because it needs to be adjusted by the operator, Impulse/P is perfectly repeatable and easy to operate. It’s also easy to clean and fast to changeover. An operator simply wipes down the stainless steel bed, snaps the decks that handle the new tablet size in place, and begins production. “We have a roller grader elsewhere in the plant. By comparison, the Impulse/P is very gentle, and it can be changed over in about half the time,” said Holwell.

Impulse/P features no sliding or rotating parts to wear. The oil-free electromagnetic drive is quiet and has no moving parts. In addition to gentle handling and high throughput, Impulse/P offers exceptional service life

and very low maintenance. “Any time we can eliminate lubricants, it’s one less thing we have to check, and it’s better for the environment,” said Holwell. “We’ve had no maintenance issues with the Impulse/P. There is nothing to maintain, really.”

With its large infeed hopper and an electronically-controlled infeed gate, Impulse/P can be easily integrated with upstream and/or downstream equipment or used as a stand-alone size grader handling batches. Jamieson uses their Impulse/P as a stand-alone unit. Producing almost 3 billion capsules and tablets per year on their 13 tablet presses that run 24 hours a day, five days a week, Jamieson’s operators assess

product coming off the presses for breakage and either run the product through Impulse/P or not, prior to tablet coating and/or packaging.

Of the 250 products that Jamieson produces, several dozen are prone to breakage and typically run through the Impulse/P. “For each product, we set up the parameters and store them in the system’s memory. Those parameters include how wide the gate opens and for how long, which regulates the product flow, because some products need to run a little slower than others,” said Bodica. Impulse/P features a touchscreen control panel with an intuitive user interface. “It’s very simple to operate.”

“In the nutraceutical industry, the quality bar is constantly rising, and we are ahead of this curve because we voluntarily follow pharmaceutical manufacturing guidelines, which are higher than nutraceutical guidelines,” noted Holwell.

“We measure the success of Impulse/P based on its ability to sort out broken tablets and fragments. Equipment uptime also matters, but product quality trumps every other consideration,” concluded Holwell. “Impulse/P does a better job removing broken tablets and fragments than manual inspection. We’ve removed the human-error factor, and done so while reducing costs.”

SYMETIX

150 Avery Street • Walla Walla, WA 99362 • USA

Phone: 800.941.3701 • Fax: 509.394.3219 • product.info@key.net • www.key.net